

## EP1

- **Cost savings—dramatically lower costs for high-throughput SNP genotyping**
- **Multiplexing—capacity for up to five reporter dyes (three standard)**
- **Flexibility—modular system design that is easily scalable**

The EP1™ system offers efficient high-throughput SNP genotyping, digital PCR and copy number variation analysis. It allows for extremely low running costs and provides the easiest workflow for low- to mid-multiplex SNP genotyping. The EP1, along with Fluidigm® integrated fluidic circuits (IFCs), streamlines the entire workflow from the setup of Dynamic Array™ IFCs and Digital Array™ IFCs to PCR thermal cycling, end-point detection and data analysis.



Designed for licensed 5' nuclease assays as well as allele-specific PCR, the EP1 allows you to easily convert from plate-based PCR to PCR with Dynamic Array IFCs without the inconvenience of switching reagents. The instrumentation consists of an end-point reader, a thermal cycler and controllers customized for your throughput requirements. As your research grows, the throughput of your EP1 system can be expanded.

Simply add additional IFC controllers and FC1™ cyclers to scale up the number of experiments per day on a single EP1. The system's economy and flexibility make it a mainstay in any research lab.

### ASSAYS

The EP1 offers an open and flexible platform to accommodate reagents and chemistries of your choice. Also, the entire system, from the footprint of the IFCs to the architecture of analysis and database software, adheres to industry standards and ensures integration with established workflows.



### ANALYSIS SOFTWARE

Bundled with data collection and analysis software for SNP genotyping and digital PCR, the EP1 SNP Genotyping Analysis software displays results in multiple formats, including scatter plots, heat maps and tabular reports. Digital PCR Analysis software allows automated analysis to easily determine copy number variations among samples.

## GENOTYPING ASSAYS

SNP Type™ assays provide a high-throughput, low-cost, single-nucleotide polymorphism (SNP) genotyping solution that enables rapid assay design and polymorphism screening. The assays are based on allele-specific PCR and combine the advantages of minimum experimental setup time and flexible assay choice with the reliability of Dynamic Array IFCs. See Figure 1.

- Designed to target species with available sequence information
- Three- to four-week design and turnaround time with customer-provided sequences (minimum of 24 assays per order)
- Access to loci-specific primer sequences assures reproducibility
- Compatible with specific target amplification (STA) protocol for improving results from samples of low quality and/or concentration, or from species with large genome sizes (>human); necessary STA primers provided

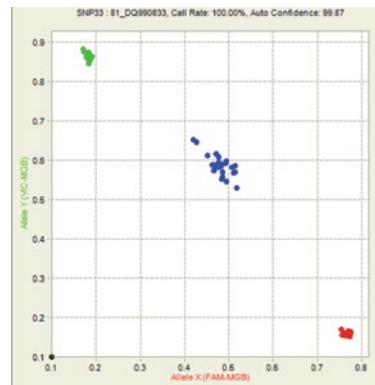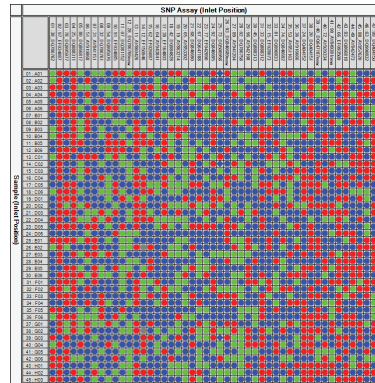

**Figure 1:**  
Call map view for 48 cattle samples and 48 SNP Type™ assays (top); cluster plot for a typical SNPtype assay (bottom)

## FC1 CYCLER

PARAMETER	
Dimensions (approx.)	9 x 8 x 19 inches (23 x 20 x 48 cm)
Software	Touchscreen interface for operation and protocol editing
Vacuum source	Internal vacuum pump
Voltage	100–230 V, 50–60 Hz
Thermal control:	
Temperature range	4–99 °C
Maximum heating rate	>5 °C/sec
Maximum cooling rate	>5 °C/sec

## IFC CONTROLLER COMPATIBILITY

PARAMETER	IFC Controller MX	IFC Controller HX	IFC Controller RX	IFC Controller WX
SNP Genotyping	48.48 Dynamic Array IFC	96.96 Dynamic Array IFC FLEXsix IFC	192.24 Dynamic Array IFC	FR48.48 Dynamic Array IFC
Digital PCR	12.765 Digital Array IFC 48.770 Digital Array IFC	—	—	—
Experiment tracking	Barcode			
Gas pressure	Internal compressor			
Interface	USB and ethernet			
IFC Controller MX, HX, RX, or WX software	Touchscreen interface for operation and sample tracking			
Dimensions (approx.)	19 x 9.5 x 13 inches; 48.5 x 24 x 33 cm			

## IFCs FOR GENOTYPING

PARAMETER	48.48 Dynamic Array IFC	96.96 Dynamic Array IFC	192.24 Dynamic Array IFC	FR48.48 Dynamic Array IFC*	FLEXsix IFC
Call rate	99.00% or greater				
Call accuracy	99.75% or greater				
Dimensions	SBS compatible (128 mm x 85 mm x 14 mm)				
Inlet spacing on input frame	4.5 mm pitch				
Liquid transfer steps	96	192	216	96	Variable
Assay inlets	48	96	24	48	6 x 12
Sample inlets	48	96	192	48	6 x 12
Reaction chambers	2,304	9,216	4,608	2,304	864
Instrument compatibility	EP1 reader, IFC Controller MX, FC1 cyclor	EP1 reader, IFC Controller HX, FC1 cyclor	EP1 reader, IFC Controller RX, FC1 cyclor	EP1 reader, IFC Controller WX, FC1 cyclor	EP1 reader, IFC Controller HX, FC1 Cyclor

\* Reusable up to five times

## IFCs FOR DIGITAL PCR

PARAMETER	12.765 Digital Array IFC	48.770 Digital Array IFC qdPCR 37k IFC
Detection sensitivity	Single copy (if copy is present in the reaction chamber)	
Dimensions	SBS-compatible (128 mm x 85 mm x 14 mm)	
Inlet spacing on input frame	4.5 mm pitch	
Minimum input volume/sample	8 µL (12 samples per array)	4 µL (48 samples per array)
Liquid transfer steps	12	48
Sample inlets	12	48
Reactions per sample	765	770
Total reaction chambers	9,180	36,960
Individual reaction volume	6 nL	0.85 nL
Total reaction volume/sample	4.6 µL (per sample)	0.65 µL (per sample)
Instrument compatibility	EP1 reader, IFC Controller MX	

## SOFTWARE SPECIFICATIONS

Fluidigm analysis software was designed to offer a simple and intuitive user interface while continuing to offer all key data analysis features for SNP genotyping and digital PCR. To simplify and expedite data analysis, the software includes these key features:

### SNP Genotyping Analysis Software

- **Automatic Cluster Calling**  
A highly accurate cluster calling algorithm automatically increases accuracy as it processes more samples.
- **Multiple Chip Scoring**  
Scoring combines multiple experiments to decrease analysis time and improve accuracy for rare genotypes.
- **Confidence Score**  
Confidence scoring automatically determines the confidence level of cluster calls.

### Digital PCR Analysis Software

- **Multiplex Capability**  
With multiplex capability, up to four reporter dyes can be used per reaction, allowing you to quantitate up to four targets per sample in each reaction.
- **Copy Number Variation (CNV)**  
CNV analysis automatically applies statistical correction and determines the ratio of your targets of interest.
- **End-Point Analysis**  
The software uses end-point image for analysis, greatly increasing the throughput of the system.

#### SYSTEM COMPONENTS

Excitation filters (center-width, in nm)	485-20, 530-20, 580-25
Emission filters (center-width, in nm)	525-25, 570-30, 630-30
Illumination	175 W Xenon Arc Lamp
Software	Fluidigm SNP Genotyping Analysis software Fluidigm Digital PCR Analysis software Fluidigm Data Collection software

#### Corporate headquarters

7000 Shoreline Court, Suite 100  
South San Francisco, CA 94080 USA  
Toll-free: 866 359 4354 | Fax: 650 871 7152  
fluidigm.com

#### Sales

North America | +1 650 266 6170 | info-us@fluidigm.com  
Europe/EMEA | +33 (0)1 60 92 42 40 | info-europe@fluidigm.com  
Japan | +81 (0)3 3662 2150 | info-japan@fluidigm.com  
China (excluding Hong Kong) | +86 (0)21 3255 8368 | info-china@fluidigm.com  
Asia | +1 650 266 6000 | info-asia@fluidigm.com  
Latin America | +1 650 266 6000 | info-latinamerica@fluidigm.com

©2014 Fluidigm Corporation. All rights reserved. Fluidigm, the Fluidigm logo, EP1, Dynamic Array, Digital Array, FC1, and SNP Type are trademarks or registered trademarks of Fluidigm Corporation in the United States and/or other countries. All other trademarks are the property of their respective owners.

**For Research Use Only. Not for use in diagnostic procedures.** PN 100-4496 B1