

Genomic DNA Preamplification Using the Advanta CFTR NGS Preamp Reagent Kit

IMPORTANT Before using this quick reference, read and understand the detailed instructions and safety guidelines in the Advanta™ CFTR NGS Library Preparation protocol for LP 48.48 IFC with Juno™ (PN 101-6270), LP 192.24 IFC with Juno (PN 101-6212), or LP 48.48 IFC with Access Array™ (PN 101-6957).

Prepare the Preamplification Reactions

IMPORTANT

- Centrifuge the enzymes before use.
- Vortex and centrifuge all other buffers and reagents before use.

- 1 In a DNA-free hood, prepare the preamplification pre-mix in a 1.5-mL tube in the order shown in Table 1.

Table 1. Preamplification pre-mix

Component		Vol. per Reaction (μL)	Vol. for 96 Reactions (μL)*
1	PCR Water (Fluidigm PN 100-5941)	1.8	216
2	4X TSP Master Mix (Fluidigm PN 101-3055)	1.25	150
3	TSP Sample Loading Reagent V2 (Fluidigm PN 101-7634)	0.25	30
4	TSP DNA Polymerase (Fluidigm PN 101-0995)	0.2	24
5	Advanta CFTR NGS Preamp Pool (Fluidigm PN 101-7279)	0.5	60
Total		4.0	480

*Includes overage.

- 2 Mix the preamplification pre-mix by briefly vortexing, and then centrifuging to bring down the contents.

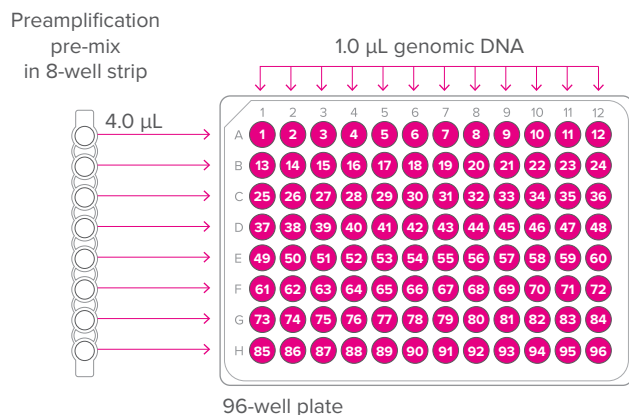


Figure 1. Preparation of preamplification reaction mix

- 3 Prepare the preamplification reaction mixes per sample as shown in Table 2 and Figure 1.

Table 2. Preamplification reaction mix

Component	Vol. per Reaction (μL)
Preamplification pre-mix (see Table 1)	4.0
Genomic DNA (gDNA), 5–30 ng/μL	1.0
Total	5.0

- a Using an 8-channel pipette, transfer preamplification pre-mix into the wells of a 96-well PCR plate.
 - b Remove the plate from the DNA-free hood and add DNA samples to each well containing pre-mix, as shown in Figure 1.
- 4 Seal the plate using an adhesive seal.
 - 5 Mix the reactions by briefly vortexing, and then centrifuge.
 - 6 Place the plate in the thermal cycler and cycle using the following thermal protocol:

Temperature	Time	Cycles	Description
95 °C	15 min		Hold
95 °C	15 sec	14	Denaturation
60 °C	8 min		Annealing/extension
4 °C	∞		Hold

Cycling time is approximately 2 hours and 45 min.

- 7 After cycling, dilute the product as shown in Table 3.

Table 3. Preamplification product dilution

Component	Vol. per Reaction (μL)
Preamplification product	2
DNA Dilution Reagent (LP 48.48: Fluidigm PN 100-9167; LP 192.24: Fluidigm PN 100-8730)	38
Total	40

- 8 Seal the plate using an adhesive seal.
- 9 Mix the diluted products by briefly vortexing, and then centrifuge.
- 10 Use the diluted preamplified gDNA as a template for preparing the sample mixes as directed in the protocol.

STOPPING POINT The diluted preamplification products can be stored at 2–8 °C if used within 1 week, or stored at –20 °C for later use.

For technical support visit fluidigm.com/support.

North America +1 650 266 6100 | Toll-free: 866 358 4354 in the US | techsupport@fluidigm.com Latin America +1 650 266 6100 | techsupportlatam@fluidigm.com
Europe/Middle East/Africa/Russia +44 1223 859941 | techsupporteurope@fluidigm.com China (excluding Hong Kong) +86 21 3255 8368 | techsupportchina@fluidigm.com
Japan +81 3 3662 2150 | techsupportjapan@fluidigm.com All other Asian countries/India/Australia +1 650 266 6100 | techsupportasia@fluidigm.com

For Research Use Only. Not for use in diagnostic procedures.

NOTICE TO PURCHASER: Limited Use License to Perform Pre-amplification with Fluidigm IFCs

The purchase of this product from Fluidigm Corporation conveys to the purchaser the limited, non-transferable right to the purchaser to use the purchased amount of this product to perform pre-amplification methods using Fluidigm instruments under license from Thermo Fisher Scientific. No right to resell this product and no other rights (such as real-time PCR methods, apparatus, reagents, or software to perform digital PCR methods) are conveyed by Thermo Fisher Scientific expressly, by implication, or by estoppel. For information on obtaining additional rights, contact outlicensing@lifetech.com or Out Licensing, Thermo Fisher Scientific, 5791 Van Allen Way, Carlsbad, California 92008.

Information in this publication is subject to change without notice. **Safety data sheet information:** fluidigm.com/sds. **Patent and license information:** fluidigm.com/legalnotices. **EU's WEEE directive information:** fluidigm.com/compliance. Fluidigm, the Fluidigm logo, Access Array, Advanta, and Juno are trademarks or registered trademarks of Fluidigm Corporation in the United States and/or other countries. All other trademarks are the sole property of their respective owners. © 2018 Fluidigm Corporation. All rights reserved. 04/2018