

Targeted DNA Sequencing Library Preparation with the LP 192.24 IFC on Juno

For safety information and detailed procedures, see the Targeted DNA Sequencing Library Preparation with Juno™ Getting Started Guide (PN 101-0414) and the Juno System User Guide (PN 100-7070).

IMPORTANT

- This document is for preparing targeted DNA sequencing libraries with Targeted DNA Seq Library Reagent kits only. To prepare sequencing libraries with Advanta™ NGS Library Prep Reagent Kits, see the Advanta NGS Library Preparation with Juno Protocol (PN 101-7878).
- Be certain that all reagents are thawed completely to room temperature and mixed thoroughly prior to use.

Obtain Target-Specific Assays

Design and order the target-specific assays ASY-MPX or ASY-MPX-P from d3.fluidigm.com. For detailed instructions about designing and ordering assays, see the D3™ Assay Design User Guide (PN 100-6812).

Dilute the Stock Assays

IMPORTANT If you are using the ASY-MPX-P diluted assay pools plate, skip this dilution step and proceed to [Prepare the Assay Pre-Mix](#).

- Immediately before use, ensure that the stock assay plate is securely sealed, and then vortex for 10–20 seconds to mix. Centrifuge the assay plate at $3,000 \times g$ for 5 minutes.
- In a DNA-free hood, dilute each pool of stock assays with DNA Suspension Buffer as calculated below. An example using the recommended volume of 2 μL for each assay from a single 96-assay pool is shown in [Figure 1](#).

Component	Starting Concentration (μM)	Number of Assays in Pool	Volume for 20X Assay Mix (μL)
Targeted DNA Seq Library Assay, one pool (Fluidigm PN ASY-MPX or ASY-MPX-P)	450/150	N	y
DNA Suspension Buffer (Teknova PN T0221)	—	—	$(180 - N) \times y$
Total	—	N	$y + [(180 - N) \times y]$

NOTE The final concentration for forward primers in each unique pool is 2.5 μM , and the final concentration for reverse primers is 833 nM.

- Using an 8-channel pipette with fresh pipette tips for each well, transfer the appropriate assay volume (y) from each column in the stock assay plate to a new 8-well strip.

- Pipet the appropriate volume of DNA Suspension Buffer into a new 1.5 mL microcentrifuge tube.
- Transfer the entire volume from each well of the 8-well strip to the tube.

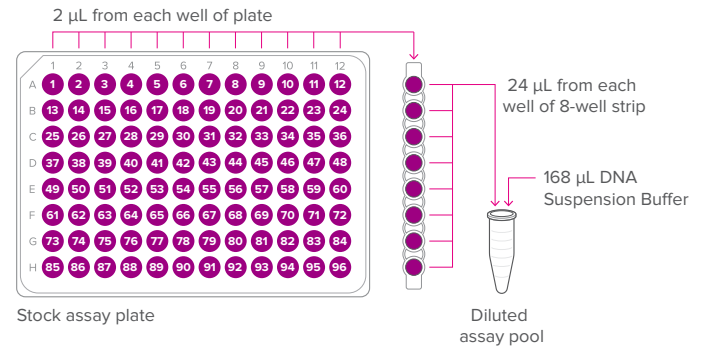


Figure 1. Example dilution of a 96-assay pool

- Repeat [step 2](#) for all other stock assay pools.
- Keep diluted assay pools at room temperature until use.

Prepare the Assay Pre-Mix

- In a DNA-free hood, combine the following components in a new 1.5 mL microcentrifuge tube:

Component	Volume per Assay Pool (μL)	Volume for 24 Assay Pools (μL)*
TSP Assay Loading Reagent (Fluidigm PN 101-0409)	2.5	75
PCR Water (Fluidigm PN 100-5941)	37.5	1,125
Total	40.0	1,200

*Includes overage. Volume is sufficient for 1 IFC. Scale up appropriately for multiple IFCs.

- Vortex the assay pre-mix for ≥ 20 seconds, and then briefly centrifuge it to bring down all components.

Prepare the 10X Assay Pools

- Vortex the diluted assay pools for 10–20 seconds to mix, and then centrifuge them at $3,000 \times g$ for 5 minutes.

- In a DNA-free hood, combine the following components in wells of a new PCR plate or in 8-well strips according to the layout shown (see Figure 2):

Component	Volume per Assay Pool (μL)*
Assay pre-mix (See Prepare the Assay Pre-Mix on page 1.)	40.0
Diluted assay pools (see Dilute the Stock Assays on page 1) or PCR water†	10.0
Total	50.0

*Includes overage.

† For unused assay inlets, replace the diluted assay pools with 10.0 μL of PCR Water.

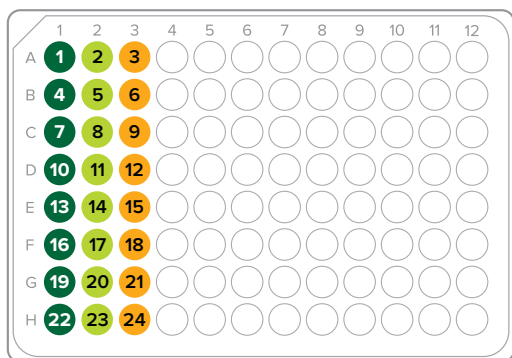


Figure 2. Layout of 10X Assay Pools

- Seal the assay plate with clear adhesive film or cap the 8-well strips.

Prepare the Sample Pre-Mix

- Manually flick the bottom of TSP DNA Polymerase tube with your forefinger to mix the enzyme (**do not vortex**). Vortex all other reagents for 20 sec, and then briefly centrifuge all reagents before use.
- In a DNA-free hood, combine in a new 1.5 mL microcentrifuge tube:

Component if using kit PN 101-5668 or 101-5669*		Volume per Reaction (μL)	Volume per IFC (μL)†
4X TSP Master Mix (Fluidigm PN 101-5786)	●	1.25	300
TSP Sample Loading Reagent (Fluidigm PN 101-0407)	●	0.25	60
TSP DNA Polymerase (PN 101-0995)	●	0.20	48
PCR Water	●	0.30	72
Total		2.00	480

*If you are using reagent kit PN 101-0406 or 101-2773, see the appendix on page 6 for the sample pre-mix volumes.

† Includes overage.

- Vortex the sample pre-mix for 10–20 seconds, and then briefly centrifuge it to bring down all components and remove bubbles.

- Pipet 56 μL of the sample pre-mix into each well of a new 8-well strip.

IMPORTANT To prevent introducing bubbles during this step, pipette only to the first stop during this transfer process. To help ensure that all liquid can be retrieved during the next step, we recommend using a microcentrifuge at maximum speed for 3 seconds.

- Keep the sample pre-mix on ice until use.

Prepare Sample Mixes

- Centrifuge the TSP Barcode Plate at 3,000 × g for 3 minutes before using.
- Obtain two 96-well plates to prepare 192 individual sample mix solutions.
- In a DNA sample hood, pipet the following components into each well of the new 96-well plates:

Component if using kit PN 101-5668 or 101-5669*	Volume per Reaction (μL)†
Sample pre-mix	2.0
Genomic DNA sample (100–200 ng/μL)	2.0
Barcode primer (Fluidigm PN 101-0744)	1.0
Total	5.0

*If you are using reagent kit PN 101-0406 or 101-2773, see the appendix on page 6 for the sample mix volumes.

† Includes overage.


- Reseal the TSP Barcode Plate. If using the barcode plate again within 2 days, store at 4 °C. Otherwise, store at –20 °C.
- Tightly seal each 96-well sample mix plate with clear adhesive film, vortex thoroughly for 20 seconds, and then centrifuge the plates at 2,500–3,000 × g for 5 minutes.
- If you observe bubbles in the wells following centrifugation, manually flick or gently snap the bottom of the affected wells with your forefinger, and then centrifuge the plates again at 2,500–3,000 × g for 5 minutes.
- Keep the sample mixes on ice until use.

Load and Run the LP 192.24 IFC

IMPORTANT

- Control line fluid on the integrated fluidic circuit (IFC) or in the inlets makes the IFC unusable. Use a new IFC.
- Vortex thoroughly and centrifuge all assay and sample mixes before pipetting into IFC inlets. Failure to do so may result in decreased data quality.
- Do not go past the first stop on the pipette.
- Before loading, ensure that there are no bubbles in the inlets.

1 Review the map for the LP 192.24 IFC:

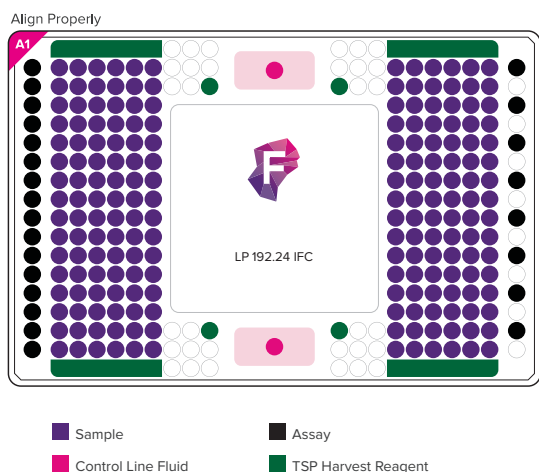

2 Pull the protective tape down and away from the bottom of the IFC. Do not turn the IFC upside down.

3 Print the life-size map of the LP 192.24 IFC, place it under the IFC, and align the notched A1 corner at the upper left of the IFC with the pink corner of the IFC map. (For a life-size map, see the Targeted DNA Sequencing Library Preparation with Juno Getting Started Guide.)


4 Load an entire syringe of Juno Control Line Fluid 192.24 into one accumulator and a second syringe into the other accumulator (pink circles on the map). Use only Juno Control Line Fluid 192.24.

5 Pipet 27 μL of TSP Harvest Reagent (Fluidigm PN 101-0743) into each of the P1, P2, P3, and P4 Inlets (green circles on the map in step 1).


6 Pipet 220 μL of TSP Harvest Reagent into each of the H1, H2, H3, and H4 harvest reservoirs (green rectangles on the map).

7 Immediately before transferring into the IFC, vortex the 10X assays and sample mixes for 10–20 seconds to mix, and then centrifuge both plates at 3,000 $\times g$ for 5 minutes. If necessary, remove any large bubbles from the wells and centrifuge again at 3,000 $\times g$ for 5 minutes.

8 Pipet 3.5 μL of each 10X assay pool into an assay inlet on the IFC (also see black circles on the map):

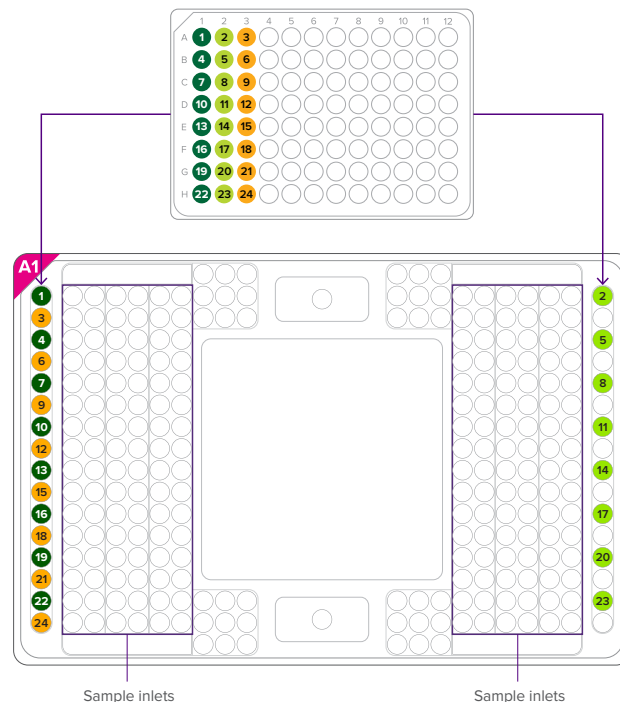

9 Pipet 3.5 μL of each sample mix into a sample inlet on the IFC (also see dark purple circles on the map). The two different colors used here illustrate the pipetting scheme:

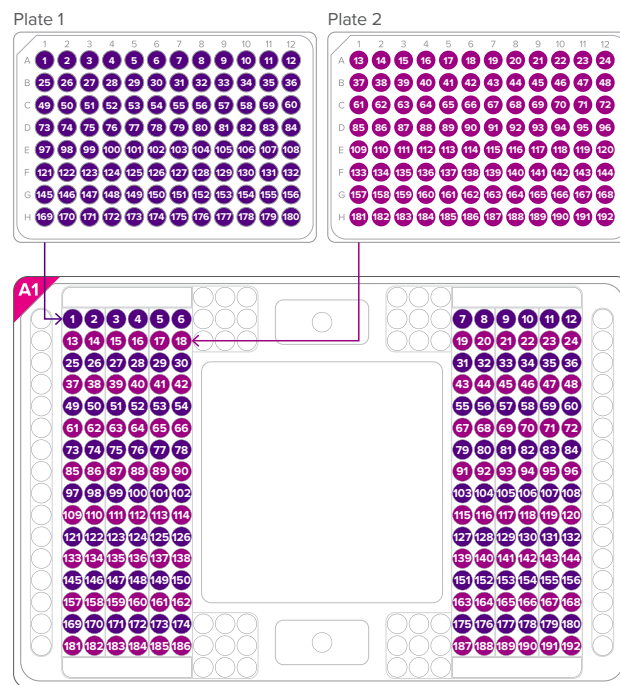

10 Cover all sample inlets with the LP 192.24 Barrier Tape with the Barrier Tape Applicator. Be sure to gently pull one of the four tabs of the barrier tape up to remove the protective sheet and expose the adhesive surface. (For detailed instructions, see the Targeted DNA Sequencing Library Preparation with Juno Getting Started Guide.)

11 Ensure that the Interface Plate TX is installed in Juno. Tap **OPEN**.

- 12** Place the IFC on the tray, and then tap **One Step LP–192.24**. Select when the script should finish (if necessary, overnight harvest delay), and then tap **RUN**.
- 13** After the IFC is finished, tap **EJECT**.
IMPORTANT Eject the IFC ≤60 minutes after the run.
- 14** After ejecting the IFC, immediately proceed to the next section.

Pool the Harvested Samples

- 1** Carefully remove the LP 192.24 Barrier Tape from the IFC by putting the IFC on a flat surface, holding the IFC with one hand, and slowly peeling off the barrier tape with the other hand
- 2** Place the life-size map of the LP 192.24 IFC under the IFC and align the notched A1 corner at the upper left of the IFC with the pink corner of the map. (For a life-size map, see the Targeted DNA Sequencing Library Preparation with Juno Getting Started Guide.)
- 3** In a post-PCR room, set an 8-channel pipette to 6.0 µL to transfer and combine the entire harvest volumes from the sample inlets of the LP 192.24 IFC directly into an 8-well strip.
IMPORTANT Be sure to transfer the entire volume from each sample inlet for best barcode uniformity of mapped reads.
- 4** Combine volumes from each 8-well strip into a single new 1.5 mL microcentrifuge tube per sample pool.
STOPPING POINT Store the 1.5 mL tube of pooled samples at 4 °C for up to one week or at –20 °C for longer storage.

First Cleanup (0.4X/0.9X Double-Sided SPRI)

- 1** Remove the Agencourt® AMPure® XP magnetic beads from 4 °C, and then leave the beads at room temperature for 30 minutes before use.
- 2** Vortex the Agencourt AMPure XP magnetic beads vigorously to ensure that they are fully suspended.
- 3** Prepare 5 mL of fresh 80% ethanol per library: Pipet 1 mL of DNase-free water into a graduated tube, and then add absolute ethanol to 5 mL. Cap the tube and mix.
- 4** Vortex the Agencourt AMPure XP magnetic beads at high speed for 1 minute.
- 5** In a new 1.5 mL microcentrifuge tube, pipet 150 µL of pooled samples. If the volume of pooled samples is <150 µL, add DNA Dilution Reagent or PCR Water to bring the volume to 150 µL.
- 6** Pipet 60 µL of AMPure XP magnetic beads into the same tube with the 150 µL of pooled samples. Expel any beads left in the pipette tip by pipetting the suspension up and down 10 times.
- 7** Vortex the suspension at high speed for 20 seconds.
- 8** Incubate the suspension at room temperature for 10 minutes.
- 9** Briefly centrifuge the tube, and then place the tube on a magnetic stand for 1–2 minutes.
- 10** Without disturbing the beads and keeping the tube on the magnetic stand, pipet the entire **supernatant** to a new tube.

- 11** Use a P10 pipette to transfer any residual volume to ensure that all supernatant has been transferred.
- 12** Pipet 75 µL of AMPure XP magnetic beads into the supernatant. Vortex the suspension at high speed for 20 seconds.
- 13** Incubate the suspension at room temperature for 10 minutes.
- 14** Briefly centrifuge the tube, and then place the tube on a magnetic stand for 1–2 minutes.
- 15** Without disturbing the beads and keeping the tube on the magnetic stand, remove and discard the supernatant.
- 16** Use a P10 pipette to remove any residual supernatant from the tube.
- 17** Keeping the tube on the magnetic stand, pipet 400 µL of 80% ethanol to wash the beads.
- 18** Incubate the tube at room temperature for 30–60 seconds.
- 19** Without disturbing the beads and keeping the tube on the magnetic stand, remove and discard the ethanol.
- 20** Repeat steps 17–19 two more times. Remove all ethanol.
- 21** Remove the tube from the magnetic stand and dry the beads: 37 °C for 1 minute, or at room temperature for 10–15 minutes.
IMPORTANT Ensure that both the beads and tube are dry before proceeding, but be careful not to over dry the beads. If over dried, beads look cracked.
- 22** To the dried beads, pipet 30 µL of DNA Dilution Reagent (Fluidigm PN 100-9167). Vortex for 20 seconds.
- 23** Incubate the suspension at room temperature for 2 minutes.
- 24** Briefly centrifuge the tube, and then place the tube on a magnetic stand for 1–2 minutes.
- 25** Keeping the tube on the magnetic stand, pipet 30 µL of the eluate to a new tube.
STOPPING POINT Store at 4 °C for up to one week or at –20 °C for longer storage.

Second Cleanup (0.8X SPRI)

- 1** Vortex the AMPure XP magnetic beads at high speed for 20 seconds.
- 2** Pipet 24 µL of Agencourt AMPure XP magnetic beads into the same tube with the 30 µL of eluate from the first cleanup. Expel any beads left in the pipette tip by pipetting the suspension up and down 10 times.
- 3** Vortex the suspension at high speed for 20 seconds.
- 4** Incubate the suspension at room temperature for 10 minutes.
- 5** Briefly centrifuge the tube, and then place the tube on a magnetic stand for 1–2 minutes.
- 6** Without disturbing the beads and keeping the tube on the magnetic stand, remove and discard the supernatant.
- 7** Use a P10 pipette to remove any residual supernatant from the tube.

- 8 Keeping the tube on the magnetic stand, pipet 190 μL of 80% ethanol to wash the beads.
- 9 Incubate the tube at room temperature for 30–60 seconds.
- 10 Without disturbing the beads and keeping the tube on the magnetic stand, remove and discard the ethanol.
- 11 Repeat steps 8–10 two more times. Remove all ethanol.
- 12 Remove the tube from the magnetic stand and dry the beads: 37 $^{\circ}\text{C}$ for 1 minute, or at room temperature for 10–15 minutes.

IMPORTANT Ensure that both the beads and tube are dry before proceeding, but be careful not to over dry the beads. If over dried, beads look cracked.

- 13 To the dried beads, pipet 30 μL of DNA Dilution Reagent. Vortex for 20 seconds.
- 14 Incubate the suspension at room temperature for 2 minutes.
- 15 Briefly centrifuge the tube, and then place the tube on a magnetic stand for 1–2 minutes.
- 16 Keeping the tube on the magnetic stand, pipet 30 μL of the eluate to a new tube.




STOPPING POINT Store the tube at 4 $^{\circ}\text{C}$ for up to 1 week or at -20°C for longer storage.

Third Cleanup (0.8X SPRI) and Quality Control

- 1 Repeat steps 1–16 in [Second Cleanup \(0.8X SPRI\)](#) with 30 μL eluate from the second cleanup. This is the purified library (before sequencing adapter is added). Store the purified library from the third cleanup at 4 $^{\circ}\text{C}$ or perform QC.
- 2 Perform QC by estimating the purified library concentration using a fluorometer and, if necessary, analyzing the purified library by Agilent® Bioanalyzer®.

Add the Sequencing Adapter to the Purified Library

- 1 In a post-PCR room, combine in a new PCR tube:

Component if using kit PN 101-5668 or 101-5669*		Volume per Reaction (μL) [†]
4X TSP Master Mix (Fluidigm PN 101-5786)		7.5
TSP Adapter Mix (Fluidigm PN 101-0408) [‡]		6.0
Purified library (before sequencing adapter is added)		4.5
PCR Water		12.0
Total		30.0

* If you are using reagent kit PN 101-0406 or 101-2773, see the appendix on [page 6](#) for the sequencing adapter volumes.

[†] Includes overage.

[‡] For dual barcoding, replace the TSP Adapter Mix with a dual index adapter mix from the Targeted DNA Seq Library Adapter Set (PN 101-2412).

- 2 Perform PCR using a stand-alone thermal cycler:

Description	Cycles	Temp.	Time
Hot start	1	95 $^{\circ}\text{C}$	15 min
PCR	10	95 $^{\circ}\text{C}$	15 sec
		60 $^{\circ}\text{C}$	90 sec
		68 $^{\circ}\text{C}$	90 sec
Final extension	1	68 $^{\circ}\text{C}$	3 min
Hold	1	4 $^{\circ}\text{C}$	∞





Clean Up the PCR Product (0.8X SPRI)

- 1 In a post-PCR room, in a new 1.5 mL microcentrifuge tube, pipet 25 μL of PCR product into 25 μL of DNase-free water. Mix, and then briefly centrifuge the tube.
 - 2 Ensure that the Agencourt AMPure XP magnetic beads are at room temperature, and then vortex the beads at high speed for 20 seconds.
 - 3 Pipet 40 μL of room temperature AMPure XP magnetic beads into the same tube with the 50 μL of diluted PCR product. (See step 1.) Expel any beads left in the pipette tip by pipetting the suspension up and down 10 times.
 - 4 Vortex the suspension at high speed for 20 seconds.
 - 5 Incubate the suspension at room temperature for 10 minutes.
 - 6 Briefly centrifuge the tube, and then place the tube on a magnetic stand for 1–2 minutes.
 - 7 Without disturbing the beads and keeping the tube on the magnetic stand, remove and discard the supernatant.
 - 8 Use a P10 pipette to remove any residual supernatant from tube.
 - 9 Keeping the tube on the magnetic stand, pipet 190 μL of 80% ethanol to wash the beads.
 - 10 Incubate the tube at room temperature for 30–60 seconds.
 - 11 Without disturbing the beads and keeping the tube on the magnetic stand, remove and discard the ethanol.
 - 12 Repeat steps 9–11 two more times. Completely remove and discard all of the 80% ethanol.
 - 13 Remove the tube from the magnetic stand and dry the beads: 37 $^{\circ}\text{C}$ for 1 minute, or at room temperature for 10–15 minutes.
- IMPORTANT** Ensure that both the beads and tube are dry before proceeding, but be careful not to over dry the beads. If over dried, beads look cracked.
- 14 To the dried beads, pipet 45 μL of DNA Dilution Reagent. Vortex for 20 seconds.
 - 15 Incubate the suspension at room temperature for 2 minutes.
 - 16 Briefly centrifuge the tube, and then place the tube on a magnetic stand for 1–2 minutes.
 - 17 Keeping the tube on the magnetic stand, pipet the entire eluate to a new, labeled tube. The eluate is the sequencing library.
- STOPPING POINT** Store the sequencing library at 4 $^{\circ}\text{C}$ for up to one week or at -20°C for longer storage.

- 18** Estimate the concentration of the sequencing library (after sequencing adapter is added) by using a fluorometer.
- 19** Compare the purified library (before sequencing adapter is added) to the sequencing library (after sequencing adapter is added) using the Agilent Bioanalyzer to ensure that the library with the sequencing adapter passes QC requirements.
- 20** Sequence the sequencing library (after sequencing adapter is added) on an Illumina® sequencer. Perform data analysis.

Appendix: Reagent Mixes for Kits PN 101-0406 and 101-2773

Sample Pre-Mix

Component		Volume per Reaction (μL)	Volume per IFC (μL)*
TSP Master Mix (Fluidigm PN 101-0994)		2.50	600
TSP Sample Loading Reagent (Fluidigm PN 101-0407)		0.25	60
TSP DNA Polymerase (PN 101-0995)		0.20	48
PCR Water		0.05	12
Total		3.00	720




*Includes overage.

Sample Mix

Component	Volume per Reaction (μL)*
Sample pre-mix	3.0
Genomic DNA sample (100–200 ng/μL)	1.0
Barcode primer (Fluidigm PN 101-0744)	1.0
Total	5.0

*Includes overage.

Sequencing Adapter

Component		Volume per Reaction (μL)*
TSP Master Mix (Fluidigm PN 101-0994)		15.0
TSP Adapter Mix (Fluidigm PN 101-0408)†		6.0
Purified library (before sequencing adapter is added)		4.5
PCR Water		4.5
Total		30.0

*Includes overage.

† For dual barcoding, replace the TSP Adapter Mix with a dual index adapter mix from the Targeted DNA Seq Library Adapter Set (PN 101-2412).

For technical support visit fluidigm.com/support.

North America +1 650 266 6100 | Toll-free (US/CAN): 866 358 4354 | techsupport@fluidigm.com
 Europe/Middle East/Africa/Russia +44 1223 598100 | techsupporteurope@fluidigm.com
 Japan +81 3 3662 2150 | techsupportjapan@fluidigm.com

Latin America +1 650 266 6100 | techsupportlatam@fluidigm.com
 China (excluding Hong Kong) +86 21 3255 8368 | techsupportchina@fluidigm.com
 All other Asian countries/India/Australia +1 650 266 6100 | techsupportasia@fluidigm.com

For Research Use Only. Not for use in diagnostic procedures.

Information in this publication is subject to change without notice. Use standard laboratory safety protocols. **Safety data sheet information:** fluidigm.com/sds. **Patent and license information:** fluidigm.com/legalnotices. **EU's WEEE directive information:** fluidigm.com/compliance. Fluidigm, the Fluidigm logo, Advanta, D3, and Juno are trademarks and/or registered trademarks of Fluidigm Corporation in the United States and/or other countries. All other trademarks are the sole property of their respective owners. © 2018 Fluidigm Corporation. All rights reserved. 06/2018